




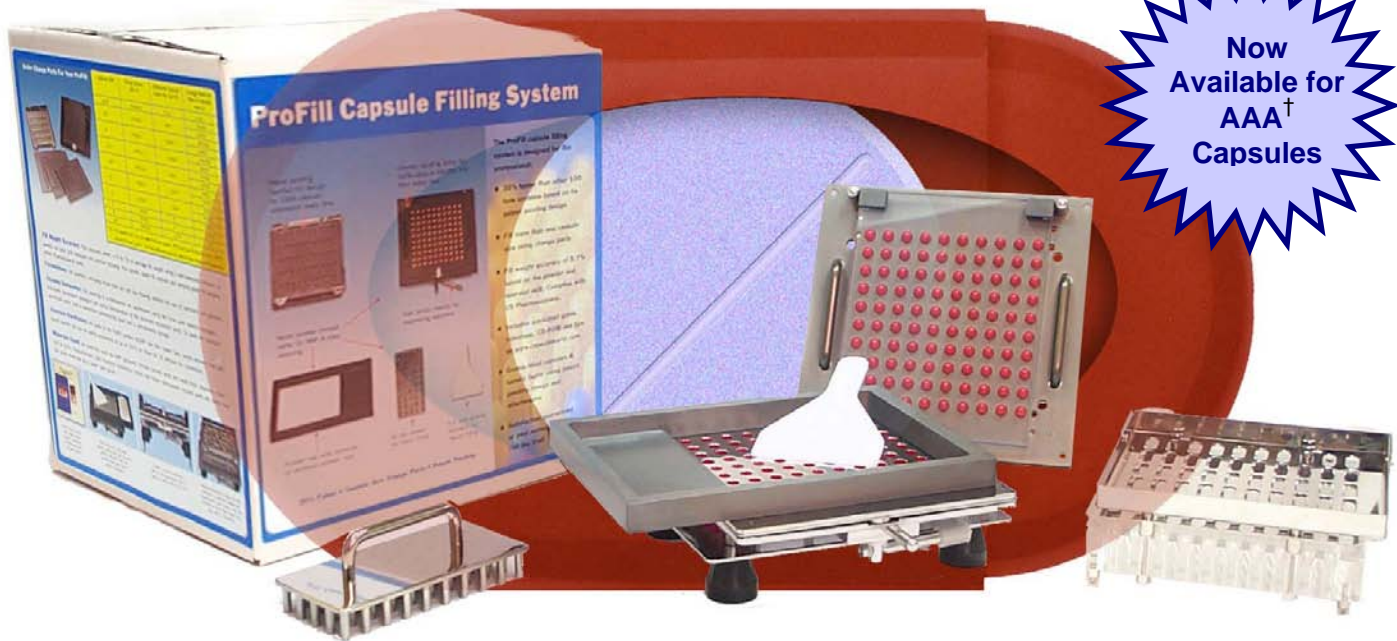

**Now Available for AAA<sup>†</sup> Capsules**





 ProFill OE (Over Encapsulator) is an alternative to hand filling small batches of DBCaps® capsules for Phase 1 & 2 studies. The Profill OE orients, separates, and locks DBCaps® capsules without vacuum or pneumatics.\*

 Double blind 1,000 to 1,500 doses/hour depending on operator speed.

 Easy to install. You can be filling DBCaps® capsules the same day you receive the ProFill OE! A great alternative to waiting for change parts or setting up ring-type or automatic fillers for a small batch.

 Available in 100 holes for sizes: AAel, AA, A, B, D; and, in 64 holes for AAA. Tablets that fit in C & E also fit in the next larger size capsule i.e. B & D.

 Special low-cost attachments make double blinding of tablets, caplets and capsules easy. For more information see reverse.

 Easy cleaning & high GMP as powder contact parts are stainless steel & anodized aluminum.

**Available for immediate delivery at US\$1850, FOB New Jersey, USA.**

\*The regular ProFill 100 mechanical separation system achieves 100% separation on standard size capsules (000-5). For the shorter DBCaps® capsules (excluding AAA) the ProFill OE incorporates a patented "single plane" capsule separation system that typically separates 95-100% of DBCaps® capsules based on capsule size and operator technique (can be as low as 90%). The remaining capsules have to be separated by hand which takes about a minute. This is preferable to separating 100% of the capsules by hand or setting up automatic filler for a small batch.

<sup>†</sup>AAA available in 64 holes. All other sizes available in 100 holes



Also available the **ProFill 100** for filling capsules from sizes 000 to 5


Visit at [www.torpac.com/100](http://www.torpac.com/100)

**Designed & Manufactured By:**

**TORPAC®**

www.torpac.com  
T: 1-973-244-1125 F: 1-973-244-1365  
333 Route 46, Fairfield, NJ 07004,



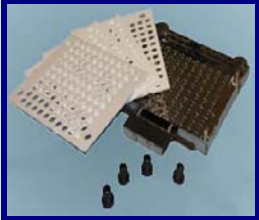

### Tablet Fill Attachment (Standard)

Over-encapsulate round tablets between 6.1-9.0mm in diameter and 2.0-5.0 mm in height. Includes stainless steel base & 4 tablet sheets for different tablets within size range given. For more information ask for the TFA Guide.

**Item #: MP TFA STD. US\$600.  
Immediate Delivery.**



### Tablet Fill Attachment (Custom)

For round tablets that are outside the size range handled by the Standard TFA. Engineering drawing of tablet required.

**Item #: MP NEC. US\$750 estimate.  
Delivery 6-8 weeks.**




### Capsule in Capsule (Standard)

Over-encapsulate filled capsules.

Capsule Size	Fits in DBcaps® size
0	AAel
1	AAel
2	AAel
3	AAel, AA, A
4	AAel, AA, A
5	AAel, AA, A, B, D

**Item #: MP NEC OCOB #\*. US\$360.  
Immediate Delivery.**  
\* # is the capsule size number

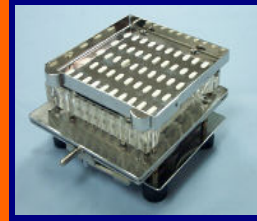

### Caplet Accessory (Custom)

For caplets or non-round tablets. Engineering drawing required.

**Item #: MP NEC. US\$750 estimate.  
Delivery 6-8 weeks.**

Attachments shown mounted on the ProFill OE Filler that is part of the ProFill OE system.

## For Filling 000-5 & el Capsule Sizes



**ProFill 100 – Hand-Held Capsule Filling System  
1500-2000 capsules/hour.**

[www.capsuleworld.com/100](http://www.capsuleworld.com/100)



**ProFill 3007 – Bench-Top Capsule Filling System  
Up to 7200 capsules/hour.**

[www.capsuleworld.com/3007](http://www.capsuleworld.com/3007)

