








-  Fast & accurate way to fill capsules without spillage.
-  Use Tamper for higher fill weights.
-  Stainless Steel (highly polished) for high GMP and easy cleanup.
-  Lightweight for use on analytical balances.
-  Available in Sizes 000 through 5, 00el, 0el; for rats 9, 9el; and, for veterinary use 11 and 12/13.



\* Free Capsules offer for sizes 000-5



**1.** Separate a single capsule. Place the capsule body (Longer, narrower part of the Capsule) in the holder.




**2.** Place the funnel over the capsule body in the holder.



**3.** Dose a pre-measured amount of powder into the capsule through the funnel, **or** put the assembly on a precision balance, tare it, and dose the powder into the capsule through the funnel.



**4.** When the required quantity of powder has been put into the funnel, use the tamper to compress filling material into capsule body.

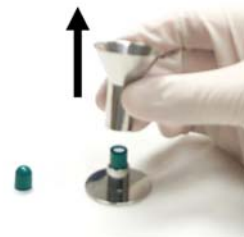

**5.** Slowly remove the funnel.



**6.** Place the capsule's cap onto the body of the capsule, pressing firmly to lock. Remove the filled capsule from the base.

### Do you need to fill more than a few capsules?

#### ProFill 100/DB



Hand-held filler up to 2000 capsules/hour. Sizes 000-5, 00el, 0el and DB.

#### ProFill 3007/8



300-hole Bench-top filler 6000 - 9000 capsules/hour. Optional vibrator & semi-automatic orienter.

#### ProCoater



Enteric / Film Coat Capsule & Caplets. Enteric coat for lower intestine drug release. Film coat for taste masking or appearance.

#### #9/9el Multi Capsule Filler



Allows you to fill 40 size 9 or 9el capsules at one time.

**TORPAC**<sup>®</sup>

www.torpac.com info@torpac.com  
T: 1-973-244-1125 F: 1-973-244-1365  
333 Route 46, Fairfield, NJ 07004, USA